


COZAP


For Coomassie Blue Removal

Destain your protein gels without changing the destain solution!

Bath Before Cozap

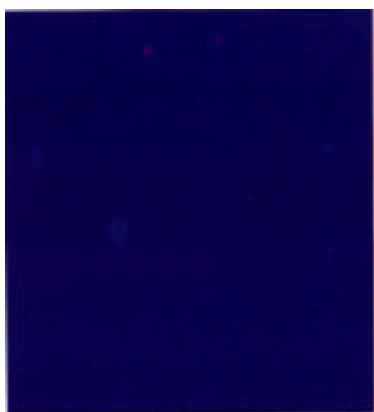

Bath After Cozap

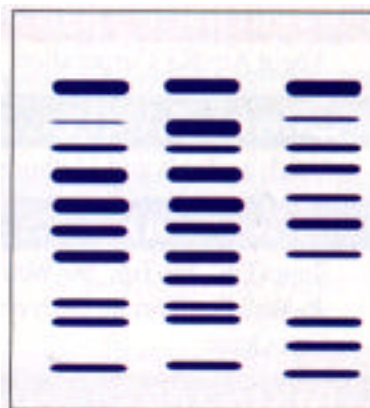

Directions:

- Place COZAP in the tank with your destaining solution. COZAP will absorb Coomassie Blue, leaving only the protein bands. (not recommended for >10% concentration of Methanol/Acetic Acid.)
- Once the gel is destained, take the gel out of the tank.
- COZAP can remain in the destaining tank until the next use.
- COZAP can be used for several gels, until it turns a deep blue color. Store COZAP at 4°C.

*For Questions or to reorder Call
The Nest Group Inc.
1-800-347-6378*

**FOR IN-VITRO USE ONLY
NOT FOR IN-VIVO, HUMAN OR ANIMAL USE**

Part #	Description	Price
SCOZ001	Pack of 25 Pads	\$85.00
SCOZ002	Pack of 100 Pads	\$307.00

- Crystal Clear Gels
- Clear Solution for Band Viewing
- Destains 5-10 mini gels each
- Faster, Cheaper, and Cleaner
- No Charcoal or Dye Residues
- No Second Wash Required
- Recovers Destaining Solution
- No Blue Bench Tops or Fingers

Cozap is a Trademark of AmiKa, Corp. patent pending.
The Nest Group, Inc. 02.03.02 S1

The Nest Group, Inc.

45 Valley Road, Southboro, MA 01772-1306

TEL#800-347-6378 or TEL# 508-481-6223

WWW URL <http://world.std.com/~nestgrp>

FAX# 508-485-5736

email: nestgrp@world.std.com