


# COZAP


## For Coomassie Blue Removal

*Destain your protein gels without changing the destain solution!*

*Bath Before Cozap*

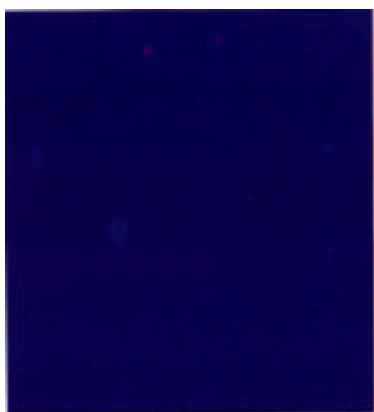

*Bath After Cozap*

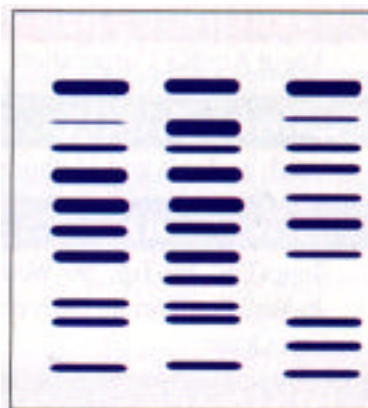

**Directions:**

- Place COZAP in the tank with your destaining solution. COZAP will absorb Coomassie Blue, leaving only the protein bands. (not recommended for >10% concentration of Methanol/Acetic Acid.)
- Once the gel is destained, take the gel out of the tank.
- COZAP can remain in the destaining tank until the next use.
- COZAP can be used for several gels, until it turns a deep blue color. Store COZAP at 4°C.

*For Questions or to reorder Call*  
The Nest Group Inc.  
1-800-347-6378

**FOR IN-VITRO USE ONLY**  
**NOT FOR IN-VIVO, HUMAN OR ANIMAL USE**

Part #	Description	Price
SCOZ001	Pack of 25 Pads	\$92.00
SCOZ002	Pack of 100 Pads	\$332.00

- Crystal Clear Gels
- Clear Solution for Band Viewing
- Destains 5-10 mini gels each
- Faster, Cheaper, and Cleaner
- No Charcoal or Dye Residues
- No Second Wash Required
- Recovers Destaining Solution
- No Blue Bench Tops or Fingers

Cozap is a Trademark of AmiKa, Corp. patent pending.  
The Nest Group, Inc. 02.03.02 S1

**The Nest Group, Inc.**

45 Valley Road, Southboro, MA 01772-1306

TEL#800-347-6378 or TEL# 508-481-6223

WWW URL <http://world.std.com/~nestgrp>

FAX# 508-485-5736

email: [nestgrp@world.std.com](mailto:nestgrp@world.std.com)